

## Environmental analysis – An electrochemical contribution




**Corrado Locati**

Electrochemistry plays a crucial role for all environment compartments, i.e., water, air and soil. In water, microbial fuel cells extract energy from wastewater, and Potentiostats monitor the associated electron transfer kinetics. In contaminated soils, Potentiostats track the redox reactions and the electrokinetic transport involved in electroremediation processes. Last but not least, Potentiostats track the electrochemical reactions that take place on catalyst surfaces – the place, where the greenhouse gas CO<sub>2</sub> is reduced to produce reactive carbon monoxide or methanol.

# Metrohm White Paper

## Water

Electrochemistry is of particular importance in characterizing microbial fuel cells, devices in which microorganisms convert oxygen and wastewater into clean water and electrical energy.

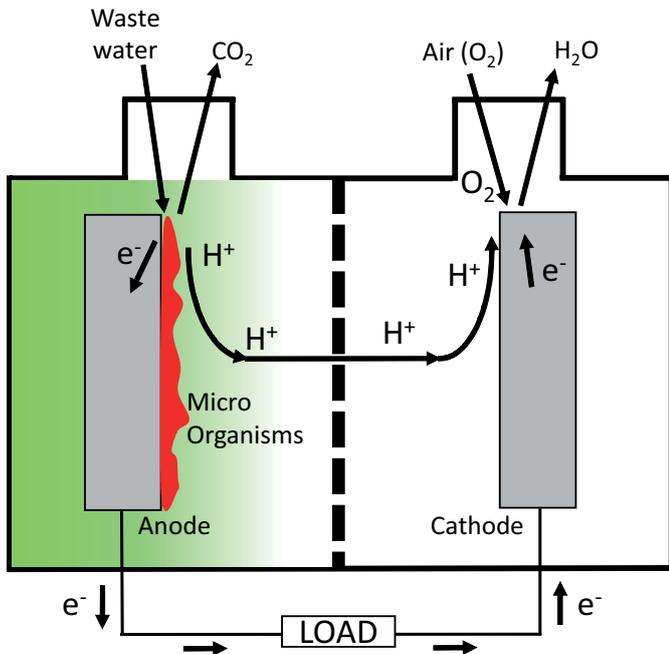


Here, the living process of microorganisms involve the breakage of complex organic molecules (the waste) into carbon dioxide ( $\text{CO}_2$ ), protons  $\text{H}^+$ , and electrons  $\text{e}^-$ . The protons cross a proton-permeable membrane (represented as a dashed vertical line), while the electrons are collected in the anode and reach the cathode through a load. At the cathode surface, the recombination of protons and electrons, together with oxygen ( $\text{O}_2$ ) from air, produces water.

In this respect, electrochemistry is an indispensable tool to provide information on the processes occurring in the microbial fuel cells [1–5]. Metrohm AUTOLAB provides complete electrochemical tools, such as the PGSTAT204 potentiostat with the FRA32M module.

## Soil

Electrochemical remediation is an in situ technique used to remove highly concentrated contaminants in soils. In a typical electrochemical remediation setup, the anode and cathode are inserted in the soil and direct current is passed through. Soil contaminants migrate at the electrodes, resulting in reduction of heavy metals ions at the cathode and oxidation of organics substances at the anode [7, 8]. The electrochemical remediation has been the first field of electrochemistry where AUTOLAB potentiostats/galvanostats have been developed. The PGSTAT302N, together with the BOOSTER20A form the best setup to perform in situ electrochemical remediation. Cyclic voltammetry, together with galvanostatic techniques, are used to monitor the redox reactions occurring in the soil under treatment.

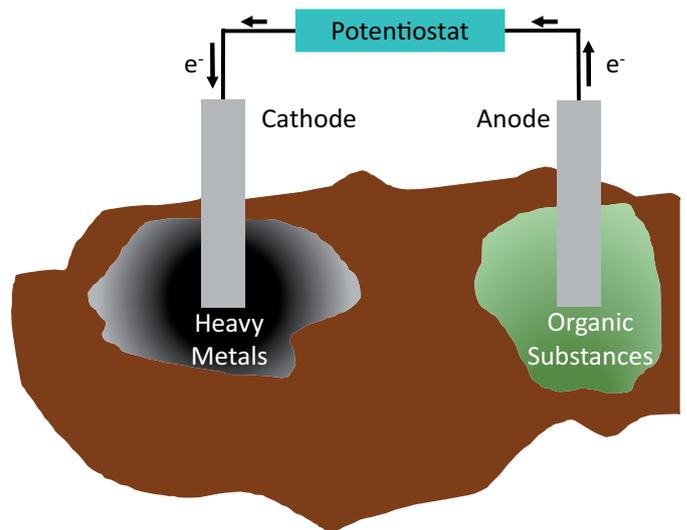


# Metrohm White Paper

## Air

One of the main pollutants affecting the atmosphere is the greenhouse gas carbon dioxide (CO<sub>2</sub>). This molecule is the main responsible for global warming, which is severely changing the climate conditions worldwide through the greenhouse effect.

Therefore, there is a need to decrease CO<sub>2</sub>, and the electrochemical reduction of carbon dioxide looks promising, since CO<sub>2</sub> is used as source for producing fuels and small organic molecules. Cyclic voltammetry or chronoamperometry are the main electrochemical techniques in use. The cathode of an electrochemical cell hosts a catalyst, which speeds up the reduction of CO<sub>2</sub>. For example, copper has been found being a good catalyst for reducing CO<sub>2</sub> dissolved in aqueous solution. In addition, other metals have been used as catalysts, as well as carbon materials [6].

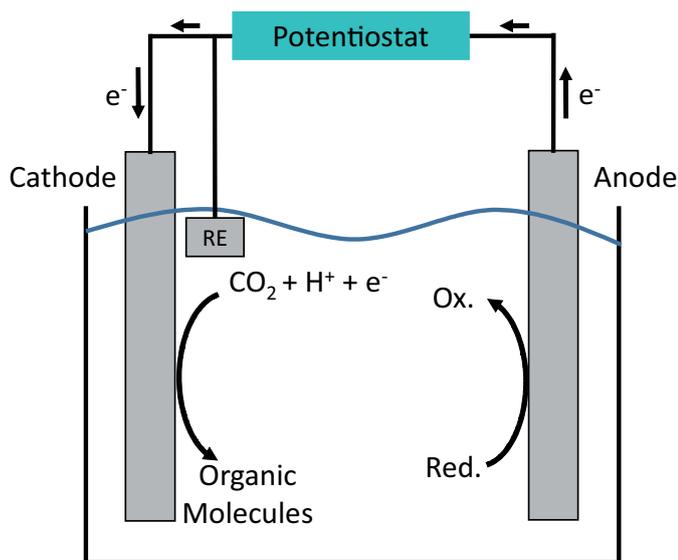

## References

- [1] B. Palys, A. Bokuna, J. Rogalski, «Poly-o-phenylenediamine as redox mediator for laccase», *Electrochimica Acta* Vol. 52 7075–7082, 2007.
- [2] R. Szamocki, V. Flexer, L. Levinb, F. Forchiasinb, E.J. Calvoa, «Oxygen cathode based on a layer-by-layer self-assembled laccase and osmium redox mediator», *Electrochimica Acta* 54 1970–1977, 2009.
- [3] C. Feng, F. Li, H. Liu, X. Lang, S. Fan, «A dual-chamber microbial fuel cell with conductive film-modified anode and cathode and its application for the neutral electro-Fenton process», *Electrochimica Acta*, Vol. 55 (6) 2048–2054, 2010.
- [4] N. Sekar and R.P. Ramasamy «Electrochemical Impedance Spectroscopy for Microbial Fuel Cell Characterization», *Microbial & Biochemical Technology*, Vol. S6 1–14, 2013.
- [5] A.K. Manohar, O. Bretschger, K.H. Nealon, F. Mansfeld, «The use of electrochemical impedance spectroscopy (EIS) in the evaluation of the electrochemical properties of a microbial fuel cell», *Bioelectrochemistry*, Vol. 72 149–154, 2008.
- [6] S. Pérez-Rodríguez, G. García, L. Calvillo, V. Celorrio, E. Pastor, and M. J. Lázaro «Carbon-Supported Fe Catalysts for CO<sub>2</sub> Electroreduction to High-Added Value Products: A DEMS Study: Effect of the Functionalization of the Support», *International Journal of Electrochemistry*, Vol. 2011 1–13, 2011.
- [7] H. Niroumand, R. Nazir, K. A. Kassim «The Performance of Electrochemical Remediation Technologies in Soil Mechanics» *International Journal of Electrochemical Science*, Vol. 7, 5708-5715, 2012.
- [8] A.T. Yeung, Y.-Y. Gu «A review on techniques to enhance electrochemical remediation of contaminated soils» *Journal of Hazardous Materials*, Vol. 195, 11–29, 2011.