


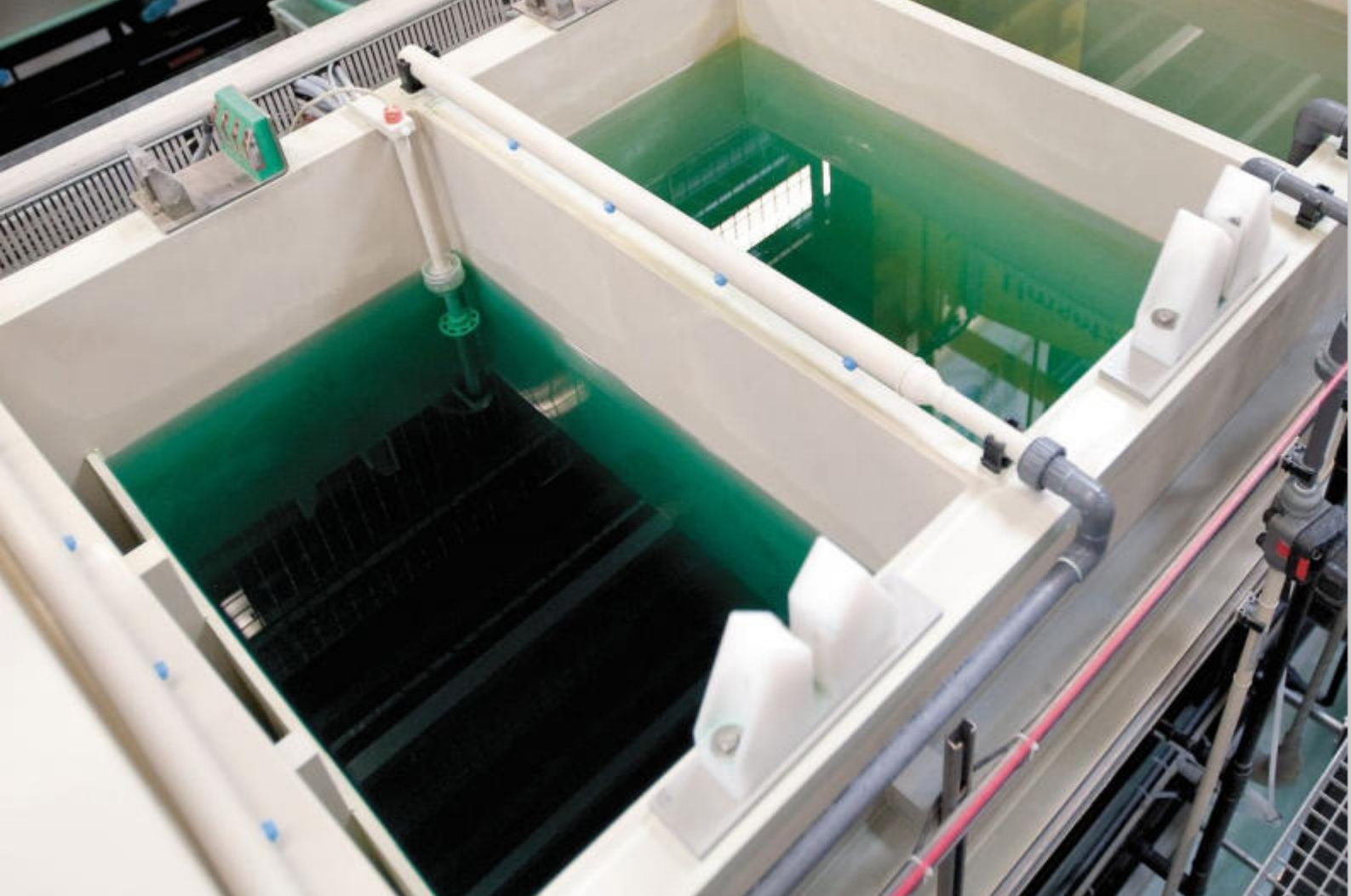

Application Note AN-V-195

無電解ニッケル浴中のよう素酸の濃度測定

Polarographic determination of iodate stabilizers

(IO_3^-)

()(PCB)ENIG()ENEPIG()



1. 884 Professional VA.

1.


Working electrode	DME
Mode	DP – Differential Pulse
Start potential	-0.1 V
End potential	-0.7 V
Peak potential iodate	-0.34 V

- :

- : Ag/AgCl/KCl (3 mol/L) . :KCl(3mol/L)

- :

mg/L

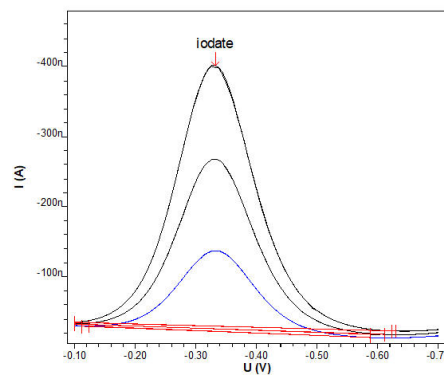

2.

2.

IO_3^- [mg/L]

2.5

Internal reference: AW CH4-0475-092007

CONTACT

143-0006 6-1-1
null 9

metrohm.jp@metrohm.jp



(MME) 884 Professional VA
semiautomated Dosinos 2

(MME) 884 Professional VA semiautomated pro
scTRACE Gold / **viva** Metrohm

CVS () CPVS () CP ()

2 800 Dosino

viva

(MME) 884 Professional VA semiautomated pro
viva



VA electrode equipment with Multi-Mode
Electrode pro for Professional VA instruments

Complete electrode set for polarographic and
voltammetric determinations. Includes Multi-Mode
Electrode pro, reference electrode, platinum
auxiliary electrode, measuring vessel, stirrer,
electrolyte solution and additional accessories for
setting up and operating the Multi-Mode Electrode.