



Application Note AN-V-230

Chromium(VI) in drinking water

Sensitive determination on the mercury film modified scTRACE Gold (DTPA method)

The guideline value for total chromium in the World Health Organization's (WHO) «Guidelines for Drinking-water Quality» is 50 $\mu\text{g/L}$. Chromium(VI) is more toxic than its trivalent form (Cr(III)) and is also less abundant. Therefore a robust and sensitive method is required to monitor its concentration in drinking water.

The mercury film modified scTRACE Gold can be used to monitor chromium(VI), offering easy handling and a high grade of stability. The sensor itself contains all necessary electrodes: the gold micro-wire working electrode, Ag/AgCl

reference, and the carbon auxiliary electrode integrated on a ceramic substrate which does not need intensive maintenance like polishing. Adsorptive stripping voltammetry (AdSV) using diethylenetriaminepentaacetic acid (DTPA) as a complexing agent and catalytic signal enhancement allows chromium(VI) detection at very low levels, with a limit of detection of 2 $\mu\text{g/L}$. The ability to replate the mercury film allows a quick and easy regeneration of the sensor. This method is best suited for field applications.

SAMPLE

Drinking water, mineral water, sea water

EXPERIMENTAL

Prior to the first determination, the ex-situ mercury film is deposited on the scTRACE Gold electrode. In the next step, the electrodes are cleaned with ultrapure water and the measuring vessel is emptied. The water sample, the supporting electrolyte with the complexing agent (DTPA) are pipetted into the measuring vessel. The determination of chromium(VI) is carried out with a 946 Portable VA Analyzer using the parameters specified in **Table 1**. The concentration is determined by two additions of a chromium(VI) standard addition solution. The scTRACE Gold is electrochemically activated prior to the first determination.



Figure 1. 946 Portable VA Analyzer

Table 1. Parameters

Parameter	Setting
Mode	DP – Differential Pulse
Start potential	-1.15 V
End potential	-1.65 V
Peak potential Cr(VI)	-1.4 V

ELECTRODES

- scTRACE Gold

RESULTS

The method is suitable for the determination of chromium(VI) concentrations up to 40 $\mu\text{g/L}$.

The limit of detection without deposition time is approximately 5 $\mu\text{g/L}$.

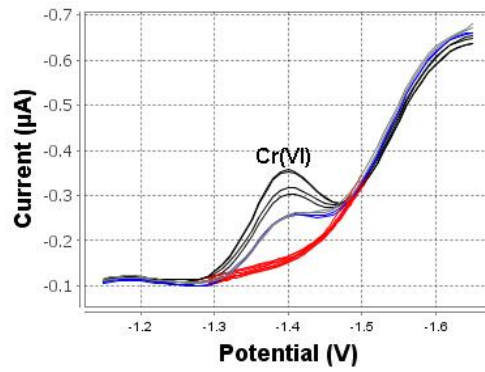


Figure 2. Determination of chromium(VI) in tap water spiked with 30 µg/L

Table 2. Result

Sample	Cr(VI) (µg/L)
Tap water spiked with 30 µg/L Cr(VI)	32.1

Internal reference: AW VA CH4-0596-042020

CONTACT

メトロームジャパン株式会社
 143-0006 東京都大田区平
 和島6-1-1
 null 東京流通センター アネ
 ックス9階

metrohm.jp@metrohm.jp

CONFIGURATION



946 Portable VA Analyzer (scTRACE Gold)

トレース範囲でヒ素、水銀、銅、鉛、亜鉛、ニッケル、コハルト、鉄、ヒスマスまたはアンチモンといった重金属測定するためのポータブル金属アナライザーです。この装置は scTRACE Gold のためのバージョンです。このシステムは、ホテニョスタットと、内蔵スターラおよび交換可能な電極かついた別個の測定スタントから構成されています。本装置は、Portable VA Analyzerソフトウェアで操作します。電源は、USBコネクタおよび内蔵の充電式電池を介して供給されます。装置は、必要なすべての付属品がキャリッジケースに入れて納品されます。