

# Differentiation of inorganic salts using Raman spectroscopy

This Application Note shows the ability of handheld Raman systems like the Mira M-1 to identify and differentiate between salts such as carbonates, phosphates, and sulfates. The main


focus of this work was to evaluate the influence of the cationic part and the water of crystallization on the Raman-spectroscopic identification of the mentioned salts.

## INTRODUCTION

Inorganic salts often have the same anionic part, while the cationic part differs. Many salts can be distinguished by the number of water molecules bond to the salt.

In this study, the influences on the Raman spectrum of the cationic part of the salt and the

number of water molecules bound to its anionic part were investigated. Although the differences between the salts are very small, the spectra that were recorded using a handheld Raman spectrometer differed sufficiently to differentiate between them.

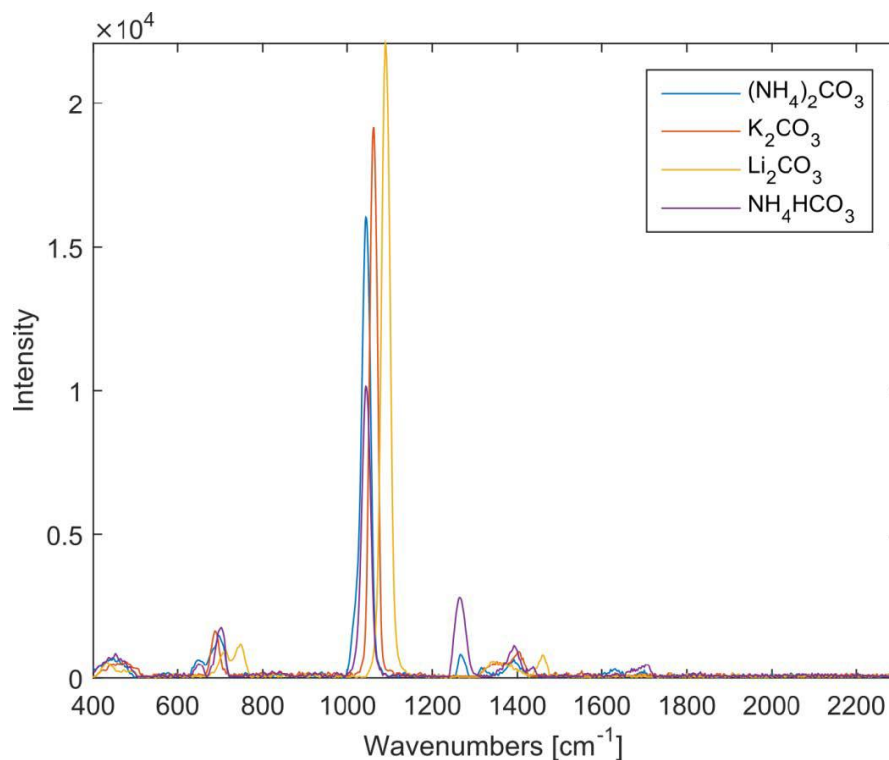

**Figure 1.** Overlaid Raman spectra of different carbonate salts.

## EXPERIMENTAL

All spectra were measured using the Mira M-1 Raman spectrometer in auto-acquisition mode, i.e., integration times were determined automatically. A laser wavelength of 785 nm and the Orbital-Raster-Scan (ORS) technique were used. The samples were analyzed in small

## RESULTS AND DISCUSSION

Figure 2 shows how the carbonate peak in the Raman spectrum shifts depending on the counter cation. This shift is significant enough to use it to differentiate between the various carbonate salts (except for ammonium carbonate and ammonium bicarbonate, where the difference is very small).

Similar results are also observed for various phosphates; the differences in the spectra are useful for their differentiation (see Figure 3) – despite the fact that there are some difficulties

vials with the vial holder adapter. Three different anions that give are Raman-active were chosen: carbonate, phosphate, and sulfate. The cationic part of the salts was varied and the resulting spectral changes were investigated.


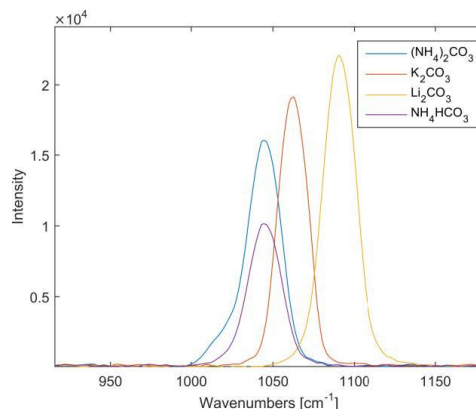


Figure 2. Main peaks of the various carbonates.


when trying to distinguish between diammonium phosphate and dipotassium phosphate (due to the similar ionic radius of potassium and ammonium).

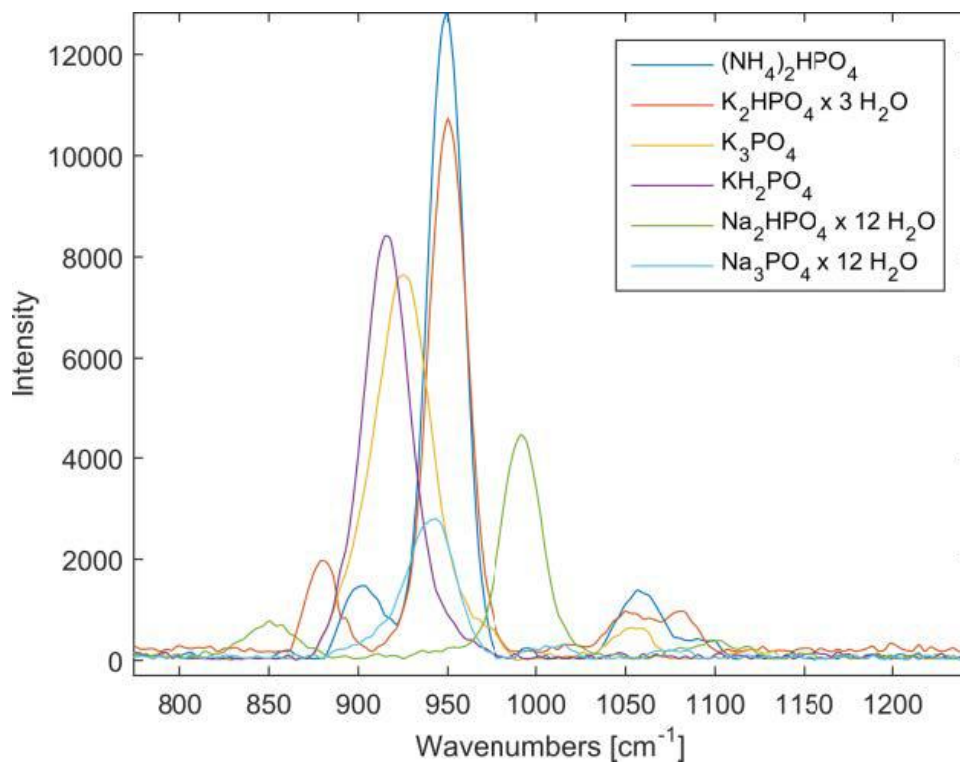

**Figure 3.** Main differences in the spectra of the phosphates.

The analyzed sulfates also differed from each other significantly, allowing their unambiguous identification with the handheld Mira M-1 (see **Figure 4**), although  $\text{CuSO}_4 \times 7 \text{H}_2\text{O}$ ,  $\text{K}_2\text{SO}_4$ ,

$\text{SnSO}_4$ , and  $\text{ZnSO}_4 \times 7 \text{H}_2\text{O}$  could not be identified unambiguously due to their similar spectra.

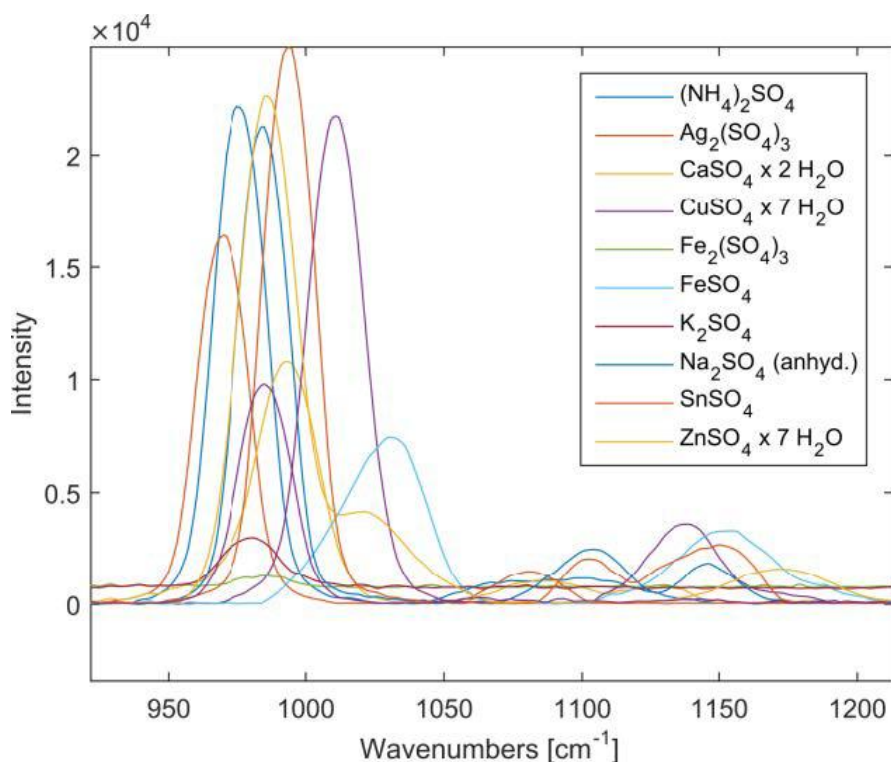

**Figure 4.** Main differences in the spectra of the sulfates.

## CONCLUSIONS

Generally speaking, handheld Raman spectrometers, such as the Mira M-1, are very useful when it comes to inorganic material identification, even when dealing with salts

having the same anionic part. Differences in the cationic parts of the salts help to unambiguously identify many salts with the Mira M-1 handheld Raman analyzer.

## CONTACT

メトロームジャパン株式会社  
 143-0006 東京都大田区平  
 和島6-1-1  
 null 東京流通センター アネ  
 ックス9階

[metrohm.jp@metrohm.jp](mailto:metrohm.jp@metrohm.jp)

## CONFIGURATION



### MIRA P Advanced

Metrohm Instant Raman Analyzer (MIRA) Pは、迅速な非破壊的計測および薬品有効成分や賦形剤などの様々な物質の検査に使用できる、高性能な携帯型ラマン分光計です。サイズはコンパクトですが、MIRA Pは非常に堅固で、弊社独自の軌道ラスタースキャン技術 (Orbital Raster Scan Technologie, ORS) を備えた作業効率の高い分光技術構造を有しています。MIRA PはFDA規則 21 CFR Part 11の要件を満たしています。

Advanced Packageには、物質を直接、またはオリジナル容器で分析することが可能なアタッチメントレンズ (レーザークラス3b)、およびガラスハイアル中のサンプル分析のためのハイアルホルターアタッチメント (レーザークラス1) が含まれています。