



Application Note AN-RS-017

Trace Detection of Thiabendazole on Bananas

Protecting consumer safety with Misa

Thiabendazole (TBZ) is a broad-spectrum pesticide used both as a fungicide in fruits and vegetables and for controlling parasites in animal feed. While some individuals experience negative side effects from consuming foods which contain TBZ, overall toxicity to humans is very low. To ensure consumer safety, regulatory agencies establish maximum residue levels (MRL) for pesticide-treated crops based on their review of risk assessment studies. For bananas, which are either aerially sprayed or dipped in

protectant solutions of TBZ, the US FDA reports a MRL of 3 $\mu\text{g/g}$, and the EU stipulates a MRL of 6 $\mu\text{g/g}$ by weight.

With Misa (Metrohm Instant SERS Analyzer), the rapid and sensitive detection of TBZ on bananas is demonstrated in formats easily adapted for food safety surveillance testing. SERS substrates for Misa offer unparalleled sampling flexibility—from dipping to swabbing, P-SERS can be adapted to the desired application.

INTRODUCTION

This application note describes the simple extraction and detection of TBZ residue on banana peels and in a processed banana

product. Samples are applied to Silver P-SERS (Ag P-SERS) strips through both **dipping** and **swab methods**.

REFERENCE SPECTRUM AND LIBRARY CREATION

To establish a reference spectrum for TBZ, a pure standard in methanol (10 μ g/mL) was analyzed using Ag P-SERS. The unique SERS spectrum

shown in **Figure 1** can be used to create a library entry for TBZ.

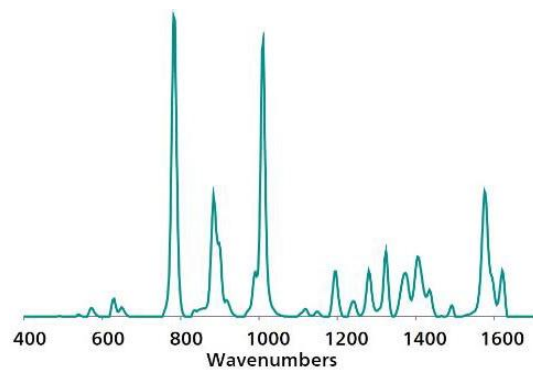


Figure 1. Standard reference Ag P-SERS spectrum of Thiabendazole.

Table 1. Experimental Parameters

Instrument		Acquisition	
Firmware	0.9.33	Laser Power	5
Software	Misa Cal V1.0.15	Int. Time	1 s
Misa Vial Attachment	6.07505.030	Averages	10
ID Kit - Ag P-SERS	6.07506.470	Raster	ON

EXPERIMENT AND RESULTS

Detection of TBZ in banana chips

Dried banana chips purchased from a local grocer were ground to a fine powder and spiked with TBZ in methanol to provide a range of test samples: 100–0.1 $\mu\text{g/g}$. Samples were mixed well and dried. Afterward, 0.1 g of each sample was added to a glass vial with 1 mL of 0.1 mol/L NaOH and shaken. After settling, an Ag P-SERS strip was **dipped** into each extract, dried for 2 minutes, and immediately placed into the Misa P-SERS attachment for measurement. **Figure 2** shows the range of detection of TBZ in ground dried banana samples from 100 $\mu\text{g/g}$ to 100 ng/g.

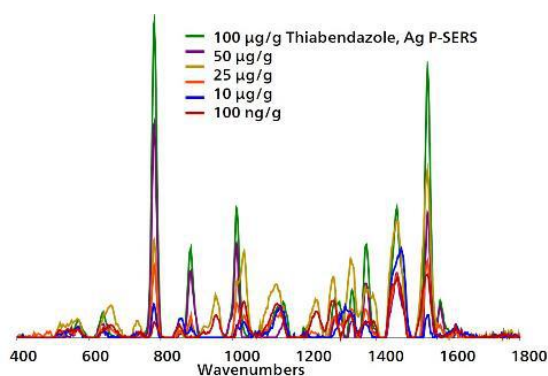


Figure 2. Overlaid baseline-corrected, dipped Ag P-SERS spectra for pulverized banana chips spiked with TBZ show detection down to 100 ng/g.

Detection of TBZ on banana peel

To simulate Ag P-SERS **swab testing**, locally purchased organic bananas were washed with tap water and isopropanol/distilled water. To simulate typical application of TBZ, a 2 cm^2 section of peel was treated with 20 μL of methanol containing different concentrations

(250, 100, 75, 50, 25, and 10 $\mu\text{g/mL}$) of TBZ. After drying, an Ag P-SERS strip wetted with methanol was used to swab the section. The strip was dried for 1 minute and placed into the Misa P-SERS attachment for measurement, with resulting spectra shown in **Figure 3**.

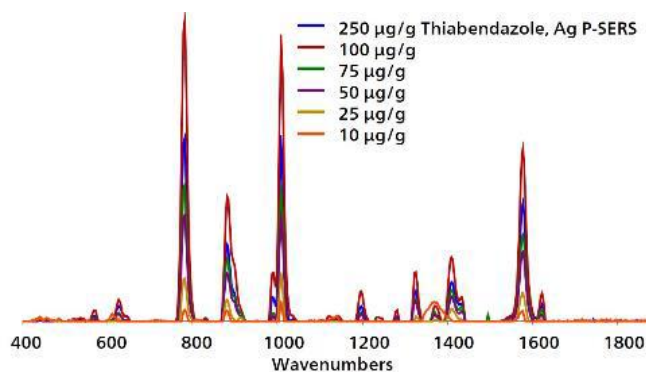


Figure 3. Overlaid baseline-corrected spectra acquired from Ag P-SERS swabs show detection of TBZ on banana peel to 100 ng/g.

Direct test of a store bought banana

In a final test, a non-organically grown unwashed banana from the grocers was **swabbed** with a methanol-moistened Ag P-SERS

strip. In **Figure 4**, detection of TBZ is confirmed by peak agreement between the spectrum from the banana swab and that of the TBZ standard.

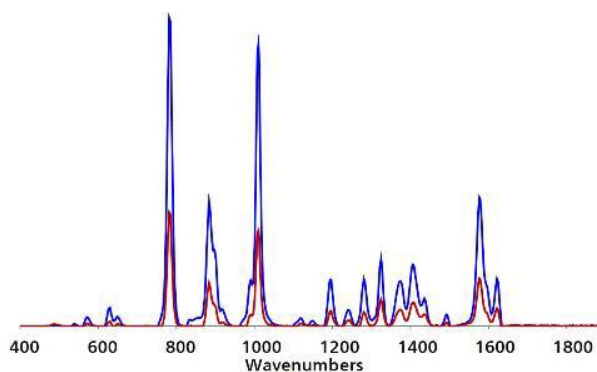


Figure 4. Direct testing of an untreated non-organic banana purchased from a grocery store returned an Ag P-SERS spectrum consistent with detection of TBZ.

FIELD TEST PROTOCOL

Detection of thiabendazole in the field

Using a disposable pipette, wet the printed end of a P-SERS strip with 2 drops of methanol. *Touch only the unprinted back side of the swab with your gloved hands.* Use the moistened, printed portion of the strip to swab the peel of a

banana. Move the strip along the peel for at least 2 cm. After drying for 2 minutes, insert the Ag P-SERS strip, with the printed portion facing down and toward the instrument, into the P-SERS attachment for measurement.

Table 2. Requirements for Field Test Protocol

ID Kit - Ag P-SERS	6.07506.470
includes:	Silver P-SERS
	Scoop
	Disposable pipettes
	2 mL glass vials
Reagents	Methanol
Test settings	Use ID Kit OP on MISA

CONCLUSION

After Misa's ability to sensitively detect TBZ on bananas is demonstrated, a successive experiment confirms post-harvest treatment of bananas with TBZ. Such fast and accurate test

results recommend Misa as an excellent solution for streamlined, onsite identification of chemically treated foods.

CONTACT

メトロームジャパン株式会社
143-0006 東京都大田区平
和島6-1-1
null 東京流通センター アネ
ックス9階

metrohm.jp@metrohm.jp

CONFIGURATION



MISA Advanced

Metrohm Instant SERS Analyzer (MISA) は、微量レベルでの違法物質、食品添加物、および食品汚染物質の迅速な検出 / 同定のための高性能な携帯可能分析システムです。MISAは、Metrohm 独自の軌道ラスタースキャン技術 (Orbital Raster Scan Technologie, ORS) を備えた高効率の分光器を有しています。これは省スペースで、より長いバッテリー寿命を持ち、現場やラボでの移動式用途にも完璧に適しています。MISA ではフレキシブルなサンプル採取を可能にする、レーザークラス1の様々なアタッチメントをご利用いただけます。アナライザーはBluetoothまたはUSBコネクタを介して操作可能です。

MISA Advanced ハッチケースは、ユーザーに Metrohmのナノ粒子溶液とP-SERSストリップを用いたSERS分析を可能にするコンフリートハッチケースです。

MISA Advanced ハッチケースには、MISAハイアルアタッチメント、P-SERSアタッチメント、ASTM校正標準、USBミニケーブル、USB電源装置、ならびにMISA装置を操作するためのMISA Calソフトウェアが含まれます。装置と付属品を安全に保管するための頑丈な保護ケースも同梱されています。



ID - P-SERS (Ag P-SERS)

IDキット - Ag P-SERS には、Mira/MisaユーザーがP-SERS銀基板でSERS分析を行うのに必要なコンポーネントが含まれています。このキットには、使い捨てのへら、滴下ヒベット、サンプルホルダー、およびP-SERS 銀ストリップが2個含まれています。