Application Note AN-T-222

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CẤU HÌNH


Eco Titrator

The compact Eco Titrator with integrated magnetic stirrer and touch-sensitive User Interface is ideal for routine analysis. It provides GLP-compliant results with minimum space requirements at all times (approx. DIN A4).

Universally compatible with almost all potentiometric titrations, such as, for example, for

- Food products: Acid content, chloride, Vitamin C, iodine and peroxide number in fats
- Water analysis: Carbonate and Ca/Mg hardness, chloride, sulfate, permanganate index
- Petrochemistry: Acid/base number, sulfide & mercaptans, chloride, bromine number
- Electroplating: Total acid, metal content, chloride
- Surfactant analysis: Anionic, cationic and non-ionic surfactants
- Photometry with the Optrode: p and m value, metals, water hardness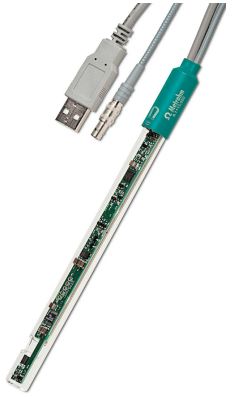

Optrode

Optical sensor for photometric titrations offering 8 different wavelengths. The wavelength can be switched using the software (tiamo 2.5 or higher) or with a magnet. The glass shaft is completely solvent-resistant and easy to clean. For example, this space-saving sensor is suitable for:

- Non-aqueous titrations in accordance with USP or EP
- Determinations of carboxyl end groups
- TAN/TBN in accordance with ASTM D974
- Sulfate determination
- Fe, Al, Ca in cement
- Water hardness
- Chondroitin sulfate in accordance with USP

The sensor is not suitable for determinations of concentrations via measurement of color intensity (colorimetry).