



Application Note AN-NIR-135

Quality control of honey with NIR spectroscopy

Simultaneous determination of color and glucose, fructose, sucrose, maltose, and turanose content with results in seconds

Honey is mainly comprised of the sugars glucose and fructose, which make up to 85% of its total weight. It additionally contains sucrose, a disaccharide composed of fructose and glucose, and other disaccharides such as maltose and turanose—present in concentrations from 0.5 to 3.5% [1]. The sugar content of honey is usually measured with high performance liquid chromatography (HPLC). Honey's color is a

quality attribute evaluated by consumers and is an important sensory property in the beekeeping market. Internationally, different types of honey are classified using the Pfund color scale. All of these honey quality parameters can be measured simultaneously in just a few seconds without any sample preparation using near-infrared spectroscopy (NIRS).

EXPERIMENTAL EQUIPMENT

Pure honey samples were measured with an OMNIS NIR Analyzer Solid (Figure 1). All measurements were performed in transflection mode (1000–2250 nm) using a 2 mm gap size reflector and 28 mm disposable vials. OMNIS Software was used for all data acquisition and prediction model development.

HPLC was the reference method used to measure the concentration of glucose, fructose, sucrose, maltose, and turanose in honey. Color was measured using a Pfund colorimeter and the Pfund scale, which ranges from 0 to 140 mm (from very light-colored honey up to the darkest honey).



Figure 1. The OMNIS NIR Analyzer Solid from Metrohm.

RESULT

The obtained NIR spectra of honey samples (Figure 2) were used to create prediction models for quantification of glucose, fructose, sucrose, maltose, and turanose content, as well as color. The quality of the prediction models was evaluated using

correlation diagrams (Figures 3–8) which display a very high correlation between the NIR prediction and the reference values. The respective figures of merit (FOM) display the expected precision of a prediction during routine analysis.

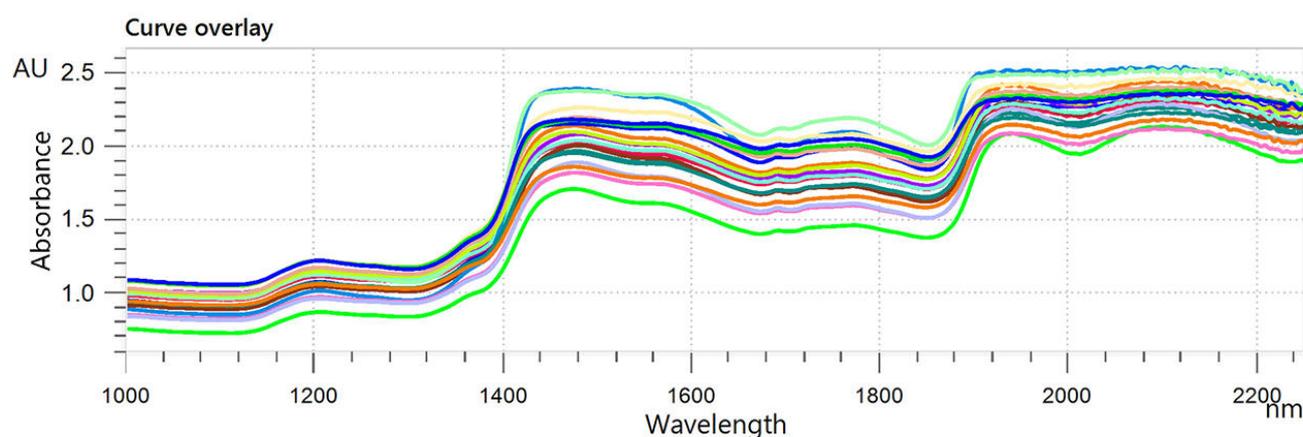


Figure 2. NIR spectra of honey analyzed on OMNIS NIR Analyzer Solid.

Result honey glucose content

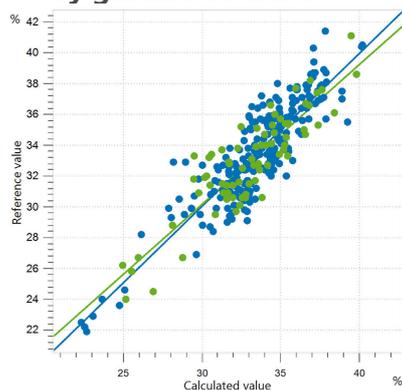


Figure 3. Correlation diagram and the respective figures of merit for the prediction of glucose content in honey. Reference values were obtained with HPLC.

R ²	SEC (%)	SECV (%)	SEP (%)
0.781	1.51	1.56	1.52

Result honey fructose content

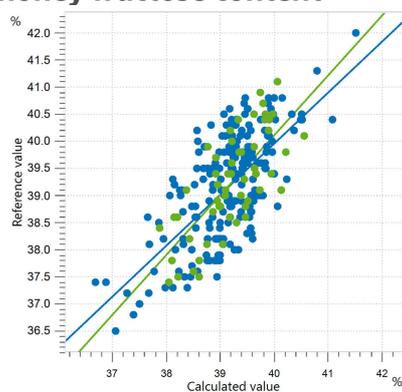


Figure 4. Correlation diagram and the respective figures of merit for the prediction of fructose content in honey. Reference values were obtained with HPLC.

R ²	SEC (%)	SECV (%)	SEP (%)
0.527	0.67	0.73	0.64

Result honey sucrose content

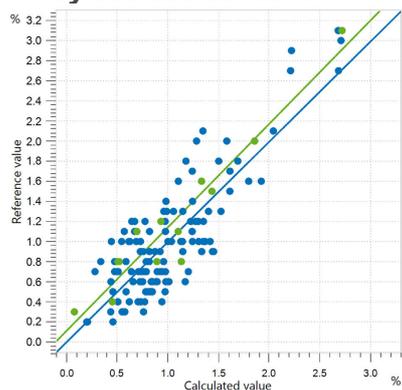


Figure 5. Correlation diagram and the respective figures of merit for the prediction of sucrose content in honey. Reference values were obtained with HPLC.

R ²	SEC (%)	SECV (%)	SEP (%)
0.917	0.29	0.32	0.25

Result honey maltose content

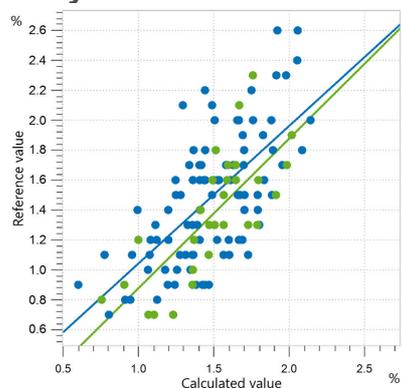


Figure 6. Correlation diagram and the respective figures of merit for the prediction of maltose content in honey. Reference values were obtained with HPLC.

R ²	SEC (%)	SECV (%)	SEP (%)
0.557	0.30	0.33	0.30

Result honey turanose content

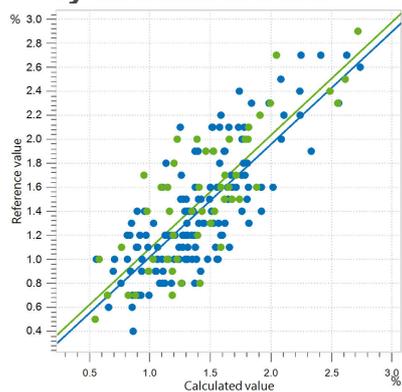


Figure 7. Correlation diagram and the respective figures of merit for the prediction of turanose content in honey. Reference values were obtained with HPLC.

R ²	SEC (%)	SECV (%)	SEP (%)
0.665	0.30	0.31	0.33

Result color

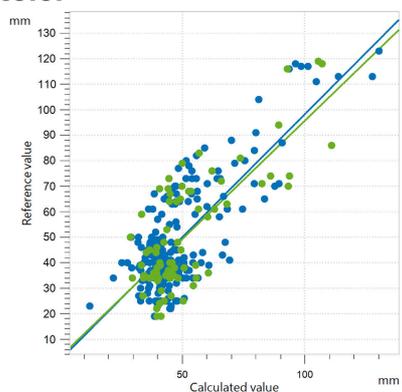


Figure 8. Correlation diagram and the respective figures of merit for the prediction of color in honey. Reference values were obtained with a Pfund colorimeter.

R ²	SEC (mm)	SECV (mm)	SEP (mm)
0.578	12.56	13.56	14.58

This Application Note displays the benefits of using near-infrared spectroscopy for quality control of honey. Color, along with glucose, fructose, sucrose, maltose, and turanose content can be measured simultaneously in only a few seconds.

Measurements performed with NIR spectroscopy

do not need any sample preparation nor solvents, saving users time and money. By using NIRS, only one analytical technology is required for sample measurement, compared to other conventional methods (**Table 1**). Finally, NIRS does not require skilled technical operators to perform the measurements, unlike HPLC.

Table 1. Overview of analytical methods used for the determination of reference values in honey.

Parameter	Method	Time to result
Glucose, fructose, sucrose, maltose, turanose	HPLC	~5 min (preparation) + ~40 min (HPLC)
Color	Pfund Method	~5 min

REFERENCE

1. Kolayli, S.; Boukraâ, L.; Sahin, H.; et al. Sugars in Honey. In *Dietary Sugars: Chemistry, Analysis, Function and Effects*; 2012; pp 3–15.

CONTACT

瑞士万通中国
北京市海淀区上地路1号院
1号楼7702
100085 北京

marketing@metrohm.com.cn

CONFIGURATION



OMNIS NIR Analyzer Solid 合固体和粘性品的近外光。

OMNIS NIR Analyzer 是一按照瑞士量准和生的近外光 (NIRS) 解决方案,用于整个生的常分析。使用新技和嵌入先 OMNIS Software 反在 NIR 光的速度、可操作性和活使用上。

OMNIS NIR Analyzer Solid 的点概:

- 在 10 秒以内量固体和粘性品
- 自化多位置量,即使在品不均匀,也能得可重的果
- 方便地嵌入自系,或者与其它分析技(滴定)
- 支持大量品容器



OMNIS NIR2 mm

液体行透射反射量的隙量 2 mm(光路度 4 mm)的反射器。

用于 28 mm 反射的一次性品瓶 (6.7402.140)。



28 mm

216 个可封的玻璃一次性品瓶,直径 28 mm,用于分析反射中的固体。用于以下分析:

- NIRS DS2500 Analyzer
- NIRS XDS RapidContent (Solid) Analyzer
- NIRS XDS MultiVial Analyzer
- NIRS XDS Masterlab Analyzer

OMNIS

A WHOLE NEW LEVEL OF PERFORMANCE

OMNIS

允机版 OMNIS 件在一台 Windows™ 计算机上行。

特性:

- 可已含有一 OMNIS 可。
- 通万通可授平台行激活。
- 不可再外计算机上使用。

OMNIS

A WHOLE NEW LEVEL OF PERFORMANCE

Quant Development

用于在独立 OMNIS Software 安装套件中写和量化模型的件可。