



Application Note AN-T-238

水中度(p和m)的定

Fast and accurate potentiometric titration of alkalinity according to ASTM D1067 and EN ISO 9963-1

概要

Alkalinity (sometimes referred to as water alkalinity) is a measure of the acid-binding capacity of water. Its degree depends on the amount of basic ions contained, mainly the carbonate content. Therefore, a distinction is made between total alkalinity (m-value) and carbonate alkalinity (p-value).

The analytical determination is carried out by titration with hydrochloric acid to pH 4.5 (turnover of the indicator methyl orange, hence «m-value»). If the initial pH of a sample is above

8.3, the acid consumption up to pH 8.3 (within the envelope of the indicator phenolphthalein, therefore «p-value») can be titrated as an intermediate step.

In this Application Note, the determination of pH and alkalinity in water with a titration method is presented. This method conforms to EPA 310.1, Standard Methods 2320 B (Titration Method), ASTM D1067, and both EN ISO 9963-1 and 9963-2.

SAMPLE AND SAMPLE PREPARATION

This application is demonstrated on tap water.

Sample preparation is not required.

EXPERIMENTAL

The determinations are carried out on an OMNIS Professional Titrator equipped with a dUnitrode with integrated Pt1000 (**Figure 1**).

An appropriate amount of water is pipetted into the titration beaker. After pH measurement, the p- and m-values are determined with fixed endpoints (FP) at pH 8.2 and 4.5 using standardized hydrochloric acid.



Figure 1. OMNIS Professional Titrator equipped with a dUnitrode with integrated Pt1000.

RESULTS

This method offers very accurate results, as displayed in **Table 1**. One exemplary curve for

alkalinity titration for water is shown in **Figure 2**.

Table 1. Results of the p alkalinity and m alkalinity of tap water expressed in mmol/L (n = 5).

Sample (n = 5)	p-value in mmol/L	m-value in mmol/L
Mean value	0.039	2.261
SD(abs)	0.002	0.001
SD(rel) in %	5.7	0.1

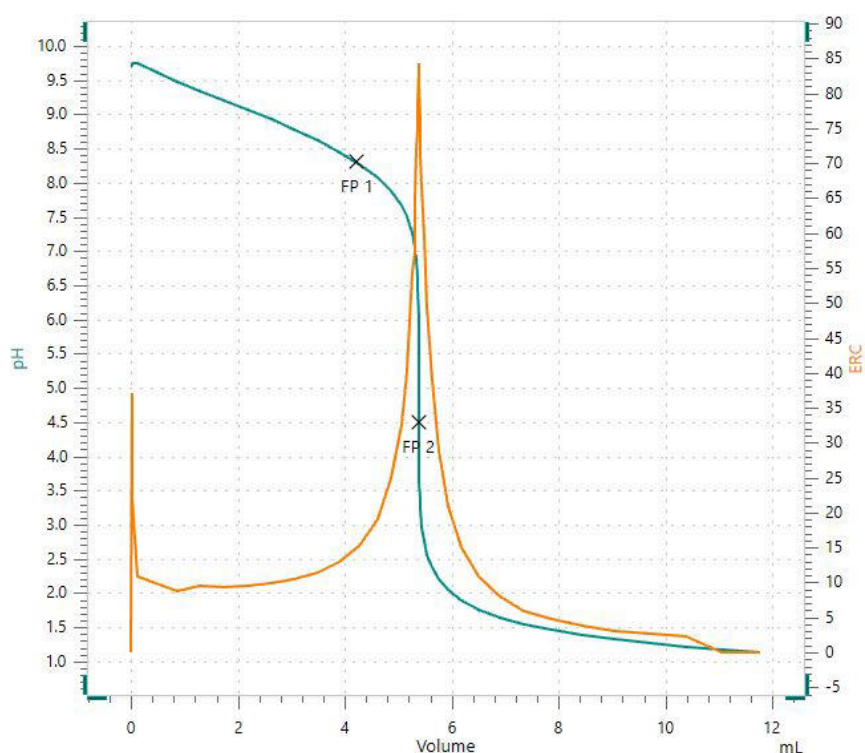


Figure 2. Titration curve showing a potentiometric determination of p alkalinity and m alkalinity with fixed endpoints at pH = 8.3 (FP1) and pH = 4.5 (FP2).

CONCLUSION

The presented titration method for alkalinity is used to accurately measure the p-value and m-value in tap water. This method conforms to EPA 310.1, Standard Methods 2320 B (Titration Method), ASTM D1067, and both EN ISO 9963-1 and 9963-2.

Precise and reliable determinations of alkalinity in water by titration with HCl are made easy using the OMNIS Professional Titrator equipped with a dUnitrode with integrated Pt1000. This system offers

users flexibility combined with high-end software. The dUnitrode is suitable for pH measurements as well as titrations in water samples. The fixed ground-joint diaphragm is resistant to contamination and the electrode works even at elevated temperatures. Aside from improving the precision and speed of the determinations, OMNIS delivers results on par with or better than other established titration systems.

CONTACT

Metrohm AG
Ionenstrasse
9100 Herisau

info@metrohm.com

CONFIGURATION



OMNIS Professional Titrator

新型、模式位分析 OMNIS Titrator,于独立行或作 OMNIS 滴定系的核心元件行,用于点和等当点滴定(一/)。由于采用 3S OMNIS Liquid Adapter 技,理化学产品从未像在一安全。可以使用量模和量管元自由配置滴定,并在需要展一台棒式拌器。包括用于使用其他滴定或加液模平行滴定的“Professional”功能可。

- 通算机或本地网控制
- 可以其他用或助溶液外接最多四个滴定模或加液模
- 棒式拌器的接方式
- 可提供不同大小的量管:5、10、20 或 50 mL
- 采用 3S 技的 OMNIS Liquid Adapter:安全理化学产品,自生厂家的原始数据

量模式和件:

- 点定滴定:“Basic” 功能可
- 点和等当点滴定(一/):“Advanced” 功能可
- 点和等当点滴定(一/),包括 5 路平行滴定:“Professional” 功能可